

Food Stamp

For food hygiene inspection and environmental surface sampling

Food Stamp RODAC (Replicate Organism Detection and Counting) plates are a *simple-to-use* bacteriological method to check CCP's (Critical Control Points) within the HACCP (Hazardous Analytical Critical Control Points) for the microbiological hygiene of surfaces.

Food Stamps are designed for hygiene control of viable bacteria on foodstuffs and working areas. It is a *simple-to-use* bacteriological method to indicate the presence of bacteria.



Easy to use

- ❖ No specific equipment required
- ❖ Easy interpretation - Specific chromogenic reaction
- ❖ Compact and easy to handle



Time saving

- ❖ No preparation needed
- ❖ Easy-to-read-results – coloured colonies





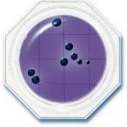
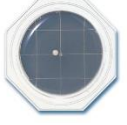
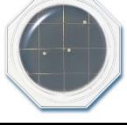

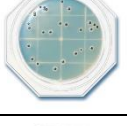
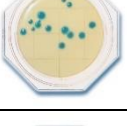
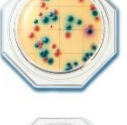
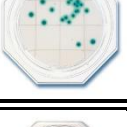
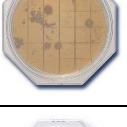

Broad applications

- ❖ Broad product menu – up to 10 different Food Stamps available for different bacteria
- ❖ Sample suitable for food stuffs, working surfaces and hygiene testing

Protocol

- Remove the cap and gently press the medium against the surface of sample. When different Food Stamps (for multiple microbes screening) are used to sample the surface, press against different areas of the sample.
- Replace the cap immediately after sampling and incubate.

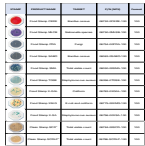

STAMP	PRODUCT NAME	TARGET	P/N (NPS)	Format
	Food Stamp CERE	Bacillus cereus	06752-0CERE-100	100
	Food Stamp MLCB	Salmonella species	06750-0MLSB-100	100
	Food Stamp PDA	Fungi	06754-00PDA-100	100
	Food Stamp SABO	Bacillus cereus	06063-0SABO-100	100
	Food Stamp SMA	Total viable count	06050-00SMA-100	100
	Food Stamp TGSE	Staphylococcus aureus	06056-0TGSE-100	100
	Food Stamp X-GAL	Coliform	06763-0XGAL-100	100
	Food Stamp XM-G	E.coli and coliform	06775-00XMG-100	100
	Food Stamp X-SA	Staphylococcus aureus	06756-00XSA-100	100
	Clean Stamp SCD*	Total viable count	06784-00SCD-100	100
	Clean Stamp SCDLP*	Total viable count	06786-SCDLP-100	100

*Clean Stamps are dedicated for Environmental Microbiology testing in Pharma setting.