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
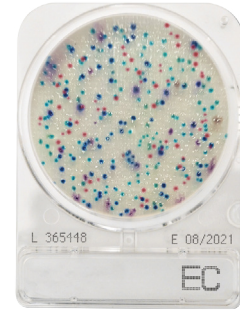
**Instructions**

Storage up to 30°C

24 months

 **Sample Preparation**

**Viable count in water or liquid foodstuff**

1 mL of sample (diluted if necessary) is deposited on the EC CompactDry plate. Pre-treat the samples according to ISO 6887 or NMKL 91.

**Viable count in solid food**

Add buffer solution (e.g. 0.9 % saline solution, PBS, buffered peptone water, etc.) to the sample and homogenize using a Stomacher® (as per manufacturer instructions). 1 mL of sample (diluted if necessary) is dispensed in the centre on the EC CompactDry plate.

**Viable count in swab test sample**

Use the swab to wipe the surface, put the swab back into the tube and screw tightly. Shake the tube to mix. 1 mL of sample (diluted if necessary) is dispensed in the centre on the EC CompactDry plate.

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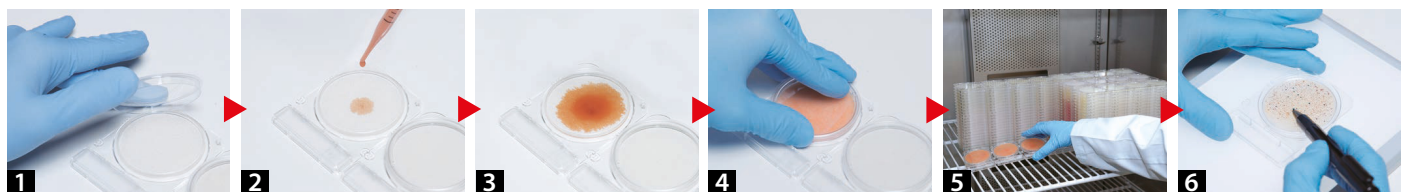
**CompactDry Protocol**

**1** Open the cap **2** Dispense 1 mL of sample on the center of the CompactDry plate OR add filter as instructed. **3** Sample diffuses automatically and evenly across the sheet transforming the dry medium into a gel within seconds. **4** Replace the cap, label and turn the plate over and **5** Incubate for the appropriate time and at required temperature. **6** Following incubation, count the number of coloured colonies.

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Incubation time 24 +/-2 hours

### Interpretation

*E.coli* forms blue colonies. Red and blue colonies together make up the total coliform group count.

### General information

*E.Coli O157* forms pink/red/purple colonies.

The growth area is 20 cm<sup>2</sup>. The back of the plate has a 1 cm<sup>2</sup> grid to ease colony counting. If there are a large number of colonies, the total viable count can be obtained by multiplying the average number of colonies in several grids by 20.

High concentrations (> 250 CFU) will cause the entire growth area to become white/pink. In this case dilute the sample.

After use please follow local disposal regulations.

### Storage and shelf life

Store at room temperature (+1 to +30 °C).

Shelf life 24 months after manufacturing.

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