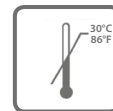
**Instructions**

Storage up to 30°C

18 months

 **Sample Preparation**

**Viable count in water or liquid foodstuff**

Pre-treat the samples according to ISO 6887 or NMKL 91.1 mL of sample (diluted if necessary) is deposited on the ETC CompactDry plate. For water samples, 100 mL or 250 mL are filtered and the membrane is incubated on CompactDry ETC.

**Viable count in solid food**

Add buffer solution to the sample and homogenize using a Stomacher® (as per manufacturer instructions). 1 mL of sample (diluted if necessary) will be dispensed in the middle on the ETC CompactDry plate.

**Viable count in swab test specimen**

Use the swab to wipe the surface, put the Use the swab to wipe the surface, put the swab back into the tube and screw tightly. Shake the tube to mix. 1 mL of sample (diluted if necessary) will be dispensed in the middle on the ETC CompactDry plate.

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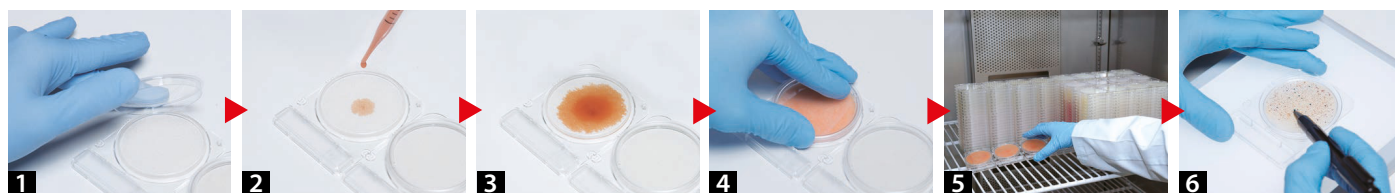
**CompactDry Protocol**

**1** Open the cap **2** Dispense 1 mL of sample on the centre of the CompactDry plate OR add filter as instructed. **3** Sample diffuses automatically and evenly across the sheet transforming the dry medium into a gel within seconds. **4** Replace the cap, label and turn the plate over **5** Incubate for the appropriate time and at required temperature **6** Following incubation, count the number of coloured colonies.

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Use the incubation time/temperature according to the national food analysis recommendations for Enterococci.

**Interpretation**

Enterococcus forms blue/blue-green colonies.

**General information**

The growth area is 20 cm<sup>2</sup>. The back of the plate has a 1 cm<sup>2</sup> grid to ease colony counting. If there are a large number of colonies, the total viable count can be obtained by multiplying the average number of colonies in several grids by 20.

High concentrations (> 250 CFU) will cause the entire growth area to become blue/blue green. In this case dilute the sample.

After use please follow local disposal regulations.

**Storage and shelf life**

Store at room temperature (+1 to +30 °C). Shelf life 18 months after manufacturing.

**Inkubationstemperatur**  
**NordVal/MicroVal: 36 +/-1°C**  
**AOAC: 37 +/-1°C**  
**Inkubationszeit 24 +/-2 Stunden**

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